


Time-of-Flight Secondary Ion Mass Spectrometry

Time-of-Flight Secondary Ion Mass Spectrometry (TOF-SIMS) is an analytical technique that uses a primary ion beam to probe the surface of a solid material.

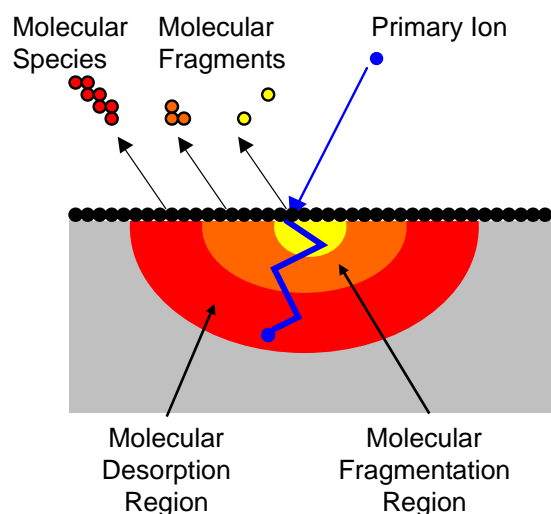

The secondary ions which desorb from the sample surface are analyzed and their mass is determined with high accuracy. As a result, elemental and molecular chemical information can be obtained on both insulating and conducting materials. TOF-SIMS allows spectroscopy for characterization of chemical composition, imaging for mapping the surface distribution of species, and depth profiling.

In the spectroscopy and imaging modes, only the outermost 1-2 monolayers of the sample surface are analyzed. To ensure the maintenance of these "static" conditions, a primary ion dose of less than 10^{12} ions/cm² is employed. Below this "static limit," roughly less than one in one-thousand surface atoms or molecules are struck by a primary ion. The actual desorption of material from the surface is caused by a "collision cascade," which is initiated by the primary ion inside the sample surface. The emitted secondary ions are extracted into the TOF analyzer by applying a potential between the sample surface and the mass analyzer. Secondary ions are generated by a pulsed primary ion source (very short pulses of <1 ns).

Secondary ions then travel through a field-free analyzer with different velocities, depending on their mass-to-charge ratio ($ke=1/2mv^2$). For each primary ion pulse, a full mass spectrum is obtained by measuring the arrival times of the secondary ions at the detector and performing a simple time-to-mass conversion. Chemical images are generated by collecting a mass spectrum at every pixel (256 x 256) as the primary ion beam is rastered across the sample surface.

TOF-SIMS is also capable of shallow depth profiling. A low-energy ion gun is operated in the DC mode for sputtering with high depth resolution. For data acquisition, the same or a different ion gun is operated in the pulsed mode. Depth profiling by TOF-SIMS allows monitoring of all species of interest simultaneously, and with high mass resolution.

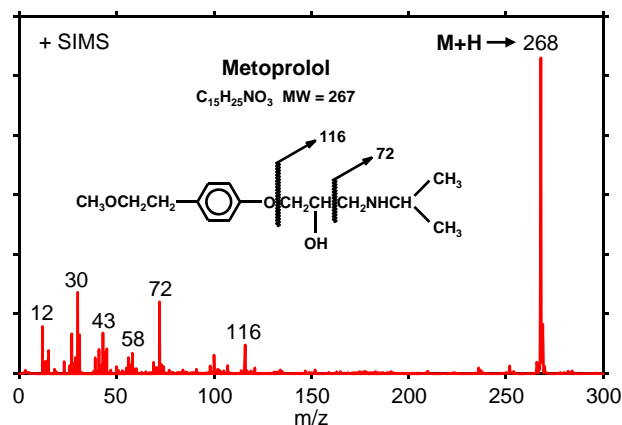


Figure 1. **Spectroscopy:** mass spectrum of Metoprolol showing molecular and fragment species.

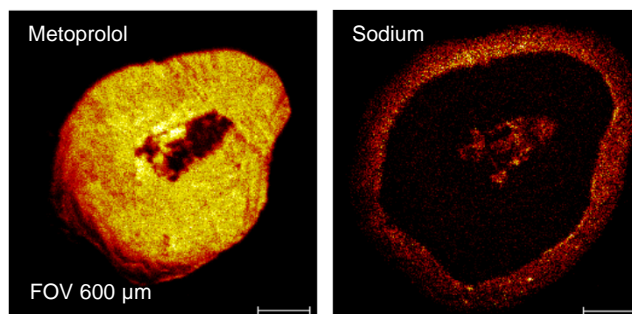


Figure 2. **Elemental and Molecular imaging:** Images of a cross-sectioned drug pellet for the molecular ion of the drug and an atomic species.

TOF-SIMS Features

- Surface Sensitivity: uppermost 20Å
- Detection limits: ppm – ppb range
- Spatial resolution: < 120 nm (determined by diameter of primary ion beam)
- Elemental and Molecular Chemical Information: practical mass range < 10,000 Da

Sample Specifications

- Sample Type: solid materials including conductors, semi-conductors and insulators
- Typical Sample Size: < 2 cm in diameter and < 10 mm thick
- Typical Analysis Area: less than 600 x 600 μm²

Typical Applications

- Polymer / Organic Coatings: on plastics, glass, metals, and paper
- Surface Contamination: additives, mold release agents, surfactants, defects
- Delamination
- Surface Modification Chemistry
- Trace Impurities
- Catalyst Surface Characterization
- Bio-materials Characterization
- Thin Film Depth Profiling

To learn more about TOF-SIMS or to discuss your specific analytical needs please contact our sales or technical staff at one of the regional support centers listed below.



Physical Electronics USA, 18725 Lake Drive East, Chanhassen, MN 55317
Telephone: 952-828-6200, Website: www.phi.com

ULVAC-PHI, 370 Enzo, Chigasaki City, Kanagawa 253-8522, Japan
Telephone 81-467-85-4220, Website: www.ulvac-phi.co.jp