



PHYSICAL ELECTRONICS

A Division of

ULVAC-PHI

SmartSoft-VP 3.2.0 for Windows 10

Improved Queue and Dashboard

SmartSoft-VP 3.2.0 is the latest release of PHI's software for VersaProbe II & III. It includes an improved queue editor for enhanced control and flexibility of data acquisition and a new sidebar dashboard supporting all options of VersaProbe III including new LEIPS, UPS, and REELS.

- Create automated analysis queues combining surveys, high-resolution scans, depth profiles, maps, and more!
- View and modify existing queue items even while the queue is running.
- Easily enable features like Z-align, pre-sputter, and use of neutralization.
- Setup advanced experiments with queue tools such as loops, pause intervals, system pressure setpoints, and more!
- Use and schedule service tasks for outgassing, startup and shutdown sequences.
- Tasks can be copied and pasted, or entire queues can be saved, loaded, and modified for routine repeated analysis.
- Queue validation feature allows user to identify and correct possible errors before starting the queue.

Queue Editor

Queue Tasks XPS Spectrum Regions

Setting
SpectrumAcq.12.26.2019.10.11.46.538 Save

| Active | Name | Sweep | Pass | Lower | Range |
|-------------------------------------|------|-------|--------|-------|-------|
| <input checked="" type="checkbox"/> | C1s | 2 | 26.000 | 280.0 | 40.0 |
| <input checked="" type="checkbox"/> | N1s | 3 | 26.000 | 396.0 | 40.0 |
| <input checked="" type="checkbox"/> | O1s | 2 | 26.000 | 528.0 | 40.0 |
| <input checked="" type="checkbox"/> | F1s | 1 | 26.000 | 684.0 | 40.0 |
| <input checked="" type="checkbox"/> | S2p | 3 | 26.000 | 165.0 | 40.0 |
| <input checked="" type="checkbox"/> | S2s | 4 | 26.000 | 225.0 | 40.0 |
| <input checked="" type="checkbox"/> | F2s | 5 | 26.000 | 25.0 | 40.0 |


Close

New Sidebar Dashboard

- Easy access to common acquisition control functions and settings
- Acquisition status can now be viewed in the dashboard to reduce the number of open windows
- Sputter Tool allows quick access to all ion gun settings
- Display of chamber pressure
- Sidebar view can be customized to minimize screen clutter

Improved Queue and Dashboard Supports all Options

- Auger Electron Spectroscopy (AES)
- Reflection Electron Energy Loss Spectroscopy (REELS)
- New automated Ultraviolet Photoelectron Spectroscopy (UPS)
- New Low Energy Inverse Photoemission Spectroscopy (LEIPS)

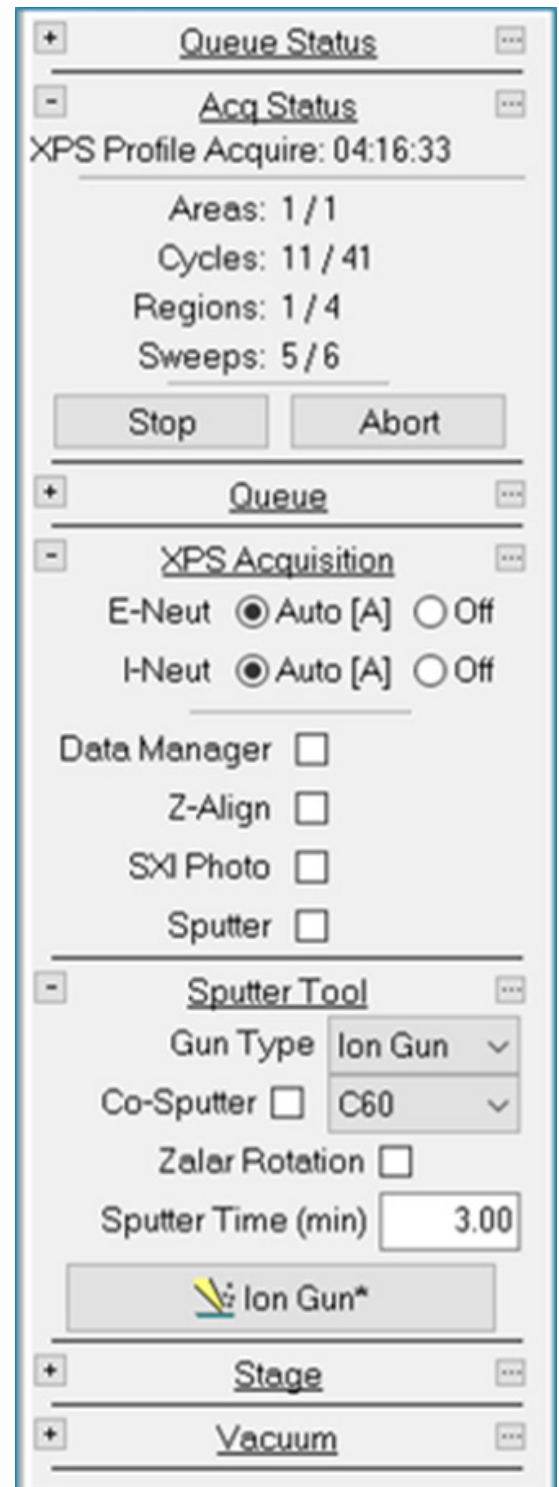

Release 3.2.0 - September 2019 (Win10: 64-bit)